


LUNA™ 3-Channel Slides

L72021 (1 box of 50 slides)
L72022 (10 boxes of 50 slides)
L72023 (10 boxes of 50 slides)
Sterile - gamma-irradiated)

Product Description

Materials	PMMA
Dimensions (W x D x H)	75 mm x 25 mm x 2.1 mm
Number of chamber	3 chambers
Chamber volume	10 μ L

LUNA™ 3-Channel Slides are individually wrapped disposable slides used with the LUNA-FX7™ Automated Cell Counter.

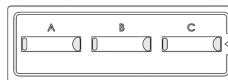

Storage

Room temperature

- Sample throughput: Up to 3 samples
- Sample loading volume: 10 μ L/chamber
- Maximum analysis volume: 1.3 μ L/chamber

Directions for Use

1. Prepare sample to count with or without staining dye.
2. Tear off the top of the wrapping and take out the slide.
3. Load 10 μ L of sample to each chamber.
4. Apply the slide to the LUNA-FX7™ Automated Cell Counter.
5. There are three-chambers to use for three sample loading.



HEADQUARTERS

FL 3
28 Simindaero 327beon-gil, Dongan-gu
Anyang-si, Gyeonggi-do 14055
South Korea

Tel: +82 (31) 478-4185

USA

7700 Little River Turnpike STE 207
Annandale, VA 22003
USA

Tel: +1 (703) 622-4660, +1 (703) 942-8867


EUROPE

1 allée Lavoisier 59650 Villeneuve d'Ascq
France

Tel: +33 (0)3 74 09 44 35

www.logosbio.com

LBSM-RD-PI-L3-001 Rev.1

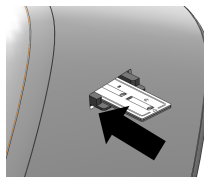

Product Care

- ! Do not press down a slide with or without wrapping.
- ! When handling, hold the edges of the wrapping, not the slide itself.
- ! Widen the opening and gently remove the slide to prevent possible scratches.
- ! Avoid fingerprints on the chamber area.
- ! Do not load any solvents buffer for counting.

Disclaimer

This product is for research use only.
Please consult the material safety data sheet for information regarding hazards and safe handling practices.

Additional information is available on our website www.logosbio.com.
© Aligned Genetics, Inc. 2021. All rights reserved.


LUNA™ 3-Channel Slides

L72021 (1 box of 50 slides)
L72022 (10 boxes of 50 slides)
L72023 (10 boxes of 50 slides)
Sterile - gamma-irradiated)

Product Description

Materials	PMMA
Dimensions (W x D x H)	75 mm x 25 mm x 2.1 mm
Number of chamber	3 chambers
Chamber volume	10 μ L

LUNA™ 3-Channel Slides are individually wrapped disposable slides used with the LUNA-FX7™ Automated Cell Counter.

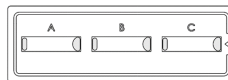

Storage

Room temperature

- Sample throughput: Up to 3 samples
- Sample loading volume: 10 μ L/chamber
- Maximum analysis volume: 1.3 μ L/chamber

Directions for Use

6. Prepare sample to count with or without staining dye.
7. Tear off the top of the wrapping and take out the slide.
8. Load 10 μ L of sample to each chamber.
9. Apply the slide to the LUNA-FX7™ Automated Cell Counter.
10. There are three-chambers to use for three sample loading.



HEADQUARTERS

FL 3
28 Simindaero 327beon-gil, Dongan-gu
Anyang-si, Gyeonggi-do 14055
South Korea

Tel: +82 (31) 478-4185

USA

7700 Little River Turnpike STE 207
Annandale, VA 22003
USA

Tel: +1 (703) 622-4660, +1 (703) 942-8867


EUROPE

1 allée Lavoisier 59650 Villeneuve d'Ascq
France

Tel: +33 (0)3 74 09 44 35

www.logosbio.com

LBSM-RD-PI-L3-001 Rev.1

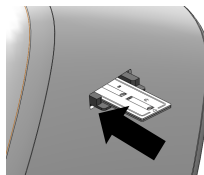

Product Care

- ! Do not press down a slide with or without wrapping.
- ! When handling, hold the edges of the wrapping, not the slide itself.
- ! Widen the opening and gently remove the slide to prevent possible scratches.
- ! Avoid fingerprints on the chamber area.
- ! Do not load any solvents buffer for counting.

Disclaimer

This product is for research use only.
Please consult the material safety data sheet for information regarding hazards and safe handling practices.

Additional information is available on our website www.logosbio.com.
© Aligned Genetics, Inc. 2021. All rights reserved.